



Instructions for LC Fittings Kit part number FT-16-8

1. Using a sharp razor or scalpel, cut the 1/16" sample tube to desired length - The cut end must be square and burr-free.
2. Slide the cap and 1/16" ferrule. Remove B.B over the cut end of the tube. Ensure that the 1/16" tube is flush or protruding slightly above the ferrule face.
3. Slide 1/8" connector over the nebulizer liquid arm, ensuring the thread is positioned as shown
4. Slip the 1/8" ferrule over the end of the nebulizer liquid arm with the tapered portion of the ferrule pointing towards the connector. Position the ferrule so that the nebulizer end protrudes slightly above the ferrule face.

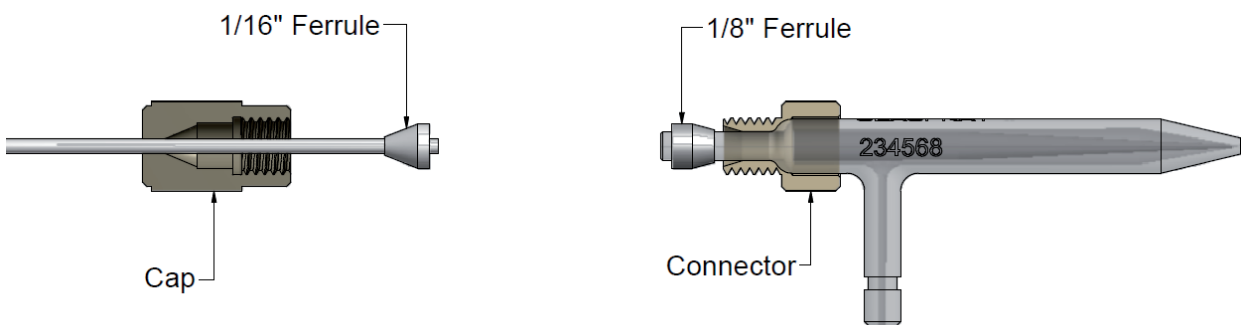

5. Hold the connector and nebulizer in one hand and start to screw together with cap until tight.
6. When finished, pull gently on the sample tube to ensure a firm connection has been made.

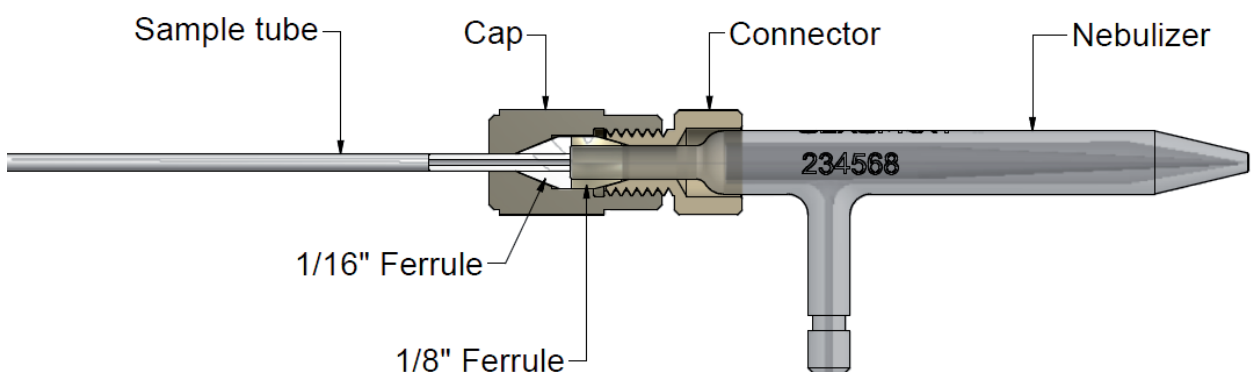

Caution! It is advised you wear safety gloves and glasses when handling glass.